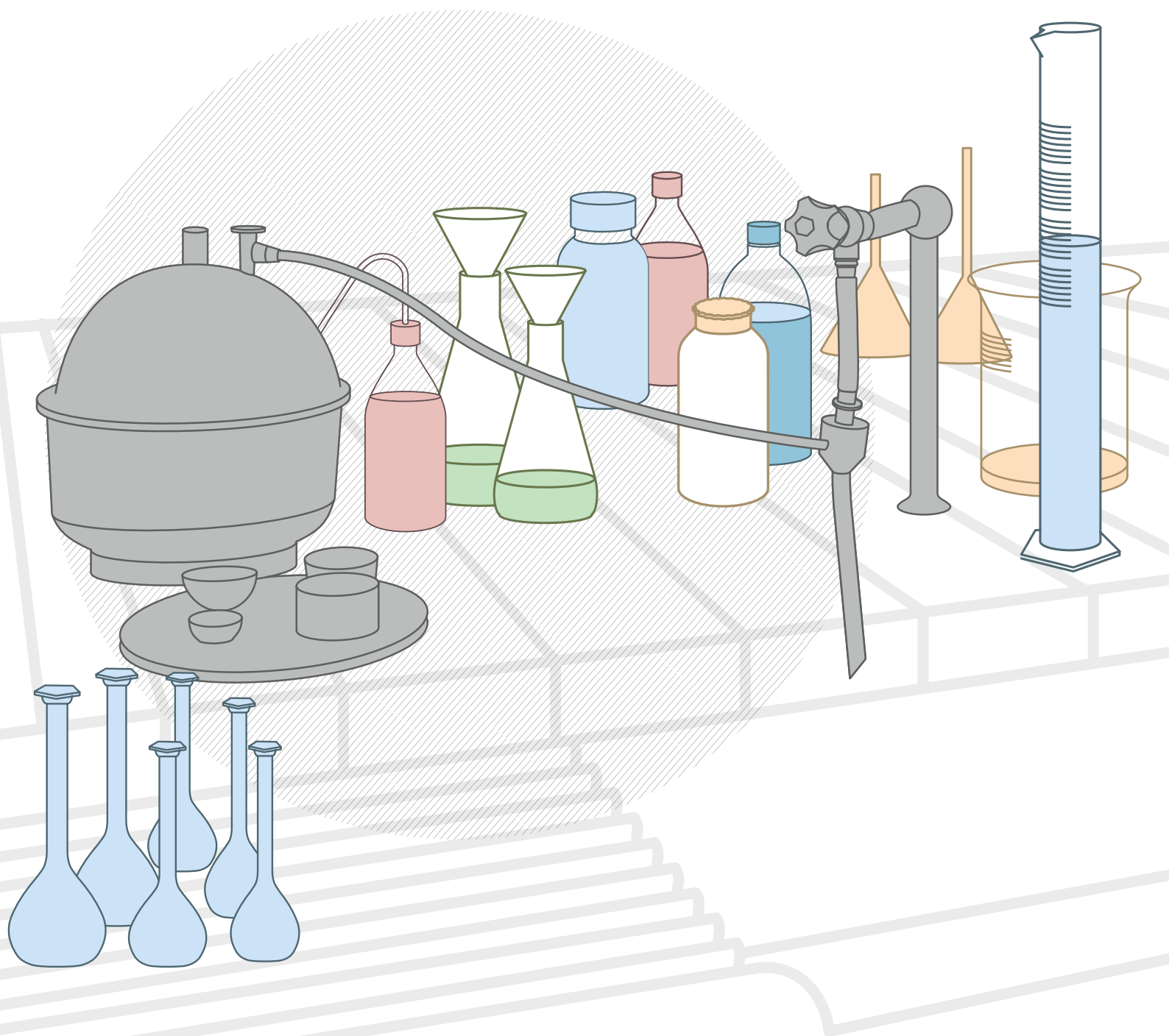
plastilab[®]

FILTRAZIONE E VUOTO - POMPE
FILTRATION AND VACUUM - PUMPS
FILTRATION ET VIDE - POMPES

**FILTRAZIONE E VUOTO
- POMPE**
**FILTRATION AND VACUUM
- PUMPS**
**FILTRATION ET VIDE
- POMPES**

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PP / PC

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**FILTRATION AND VACUUM
DESICCATORS**

PP / PC

Implosion proof desiccators. Will take vacuum of -740 mm Hg with loss (absolute zero: -760 mm/HG) not exceeding 20 mm Hg in 24 hours. The transparent polycarbonate cover seals onto a neoprene 'O' ring set into the flange of the PP base. A little grease spread on the 'O' ring will assist in establishing initial vacuum. Within the desiccator is a removable pan for holding a drying agent. Vacuum release tap/non return valve assembly is included; specially grooved stopper in this assembly allows air, when required, to be re-admitted very slowly to avoid disturbing the contents.

Usage Instructions:

- remove the cover in PC - insert the samples in the pan
- open the vacuum retention valve (PE needle) connect the tube of the pump to the desiccator spigot in order to create vacuum close the vacuum retention valve

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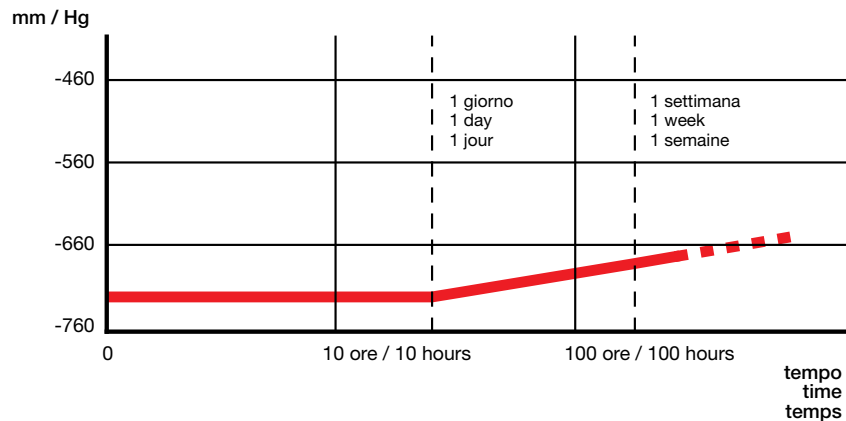
DESICCATOR PLATES

PP

These plates made of PP are extremely useful as a support for crucibles, Petri dishes etc. Use at room temperature. Not suitable for use with hot crucible.

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**SPARE VACUUM RETENTION VALVE
PC WITH PE NEEDLE**

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
This pump will work on pressures up to 10 Kg/cm². Built in non return valve to eliminate back flow. Useful for sucking to waste any split chemicals as well as for filter work. Easily dismantled for cleaning. Autoclavable.

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The graph below details the performance of the pump under different conditions. Example: Operating pressure of water: 1 Kg/cm². Attained Vacuum: 15 mm/Hg. Will evacuate a 5 litre container in about 20 minutes.

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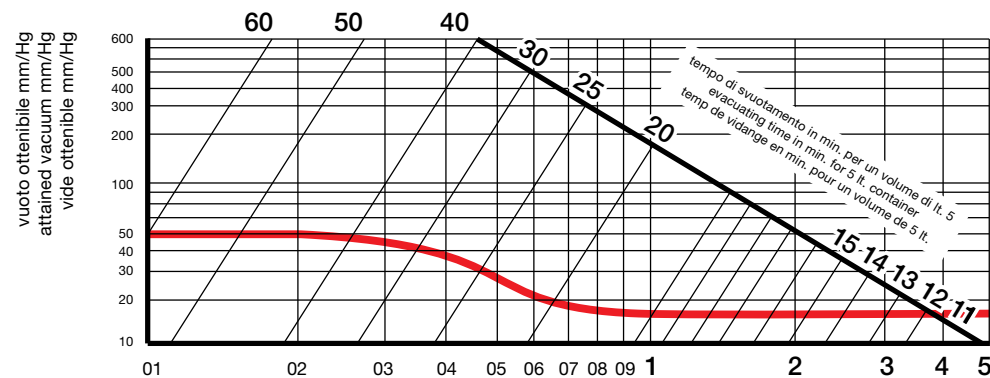NON-RETURN VALVE

PP INNER DISK: DUTRAL

Prevents water back-flow when pressure falls. Useful for any tubing system and vacuum line. Maximum safe working pressure: 2 BAR (equal approx. 28.4 P.S.I.). Minimum safe working pressure 0,07 BAR at room temperature (20°C).

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FUNNELS, BUCHNER

PP


Robust, lightweight. Unbreakable and resistant to corrosion and heat. Two piece construction. Top and bottom can be separated for easy and total cleaning. Autoclavable. Excellent chemical resistance. **SUITABLE FOR FOODSTUFF.**

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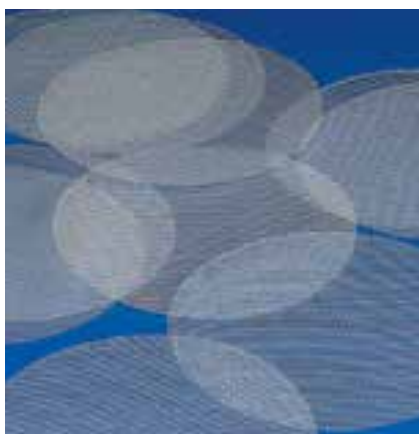DISKS FOR BUCHNER FUNNELS

PE

Useful interface between filter paper and perforated Buchner plate. Avoids clogging of filter holes. Texture 0,5 mm.

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HAND OPERATED VACUUM PUMP

HIGH IMPACT PS

Light, portable hand operated vacuum pump. Attains a vacuum of 625 mm Hg (25"). Pumping rate is 15 ml per stroke. Built in vacuum release valve. Can be used as pressure pump by transferring release valve cap to vacuumport. Nozzle fits standard 6 mm (1/4") i.d. tubing.

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ANALYTICAL FUNNELS

PP

Designed specifically for analytical chemistry. Each size takes a standard size filter paper. Body of funnel is exact 60° angle. Provided with external and internal ribs. Autoclavable. **SUITABLE FOR FOODSTUFF.**

Art.	Ø mm	Cap. ml	Lung. gambo mm Stem length mm Long. tige mm	Gambo Ø mm Stem Ø mm Tige Ø mm	Tot. L mm
152	100	200	82	11	161,5
153	120	350	86	11	184
155	150	700	115	14	238
156	183	1250	140	14	293

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LONG-STEM FUNNELS

PP

The 60° angles sides are provided with ribs for fast filtration and allow the use of standard size filter papers. **SUITABLE FOR FOODSTUFF.**

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VISCOUS LIQUID PUMPS

PP

Dispensing pump made of PP operated by a unique lever-action stroke. Delivery of approximately 10 ounces (ab. 300 ml) per stroke, with capability of 6 gallons per minute (ab. 22-23 litres/min). Will handle the most viscous materials (equivalent to SAE 90 oil). Spigot orientation of 360° allows versatility in dispensing. Complete with shank.

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